Originalbetriebsanleitung
Translation of the original Operating Instruction

Steripur 417

2/2 Wege-Membranventil

2/2 Way Diaphragm Valve

Inhalt

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Content

1.	Introduction.....	24
1.1.	Information about the operating instructions.....	25
2.	Safety.....	26
2.1.	Safety Information.....	26
2.2.	Hazard classification	26
2.3.	Intended use.....	27
2.4.	Misuse	27
2.5.	General safety provisions	27
2.6.	Residual risks	28
2.7.	Responsibilities of the operator	28
2.7.1	Appointment and instruction of responsible persons.....	28
2.7.2.	Information obligation	28
2.8.	Description of target groups.....	29
3.	Transport / Storage / Disposal	29
3.1.	Delivery.....	29
3.1.1.	Scope of delivery.....	29
3.1.2.	Notification about damage	29
3.2.	Transport	29
3.3.	Storage.....	30
3.4.	Disposal.....	30
4.	Technical Data	31
4.1.	General technical data.....	31
4.2.	Operating conditions	32
4.2.1.	Operating temperatures	32
4.2.2.	Operating pressure	32
4.2.3.	Control pressure	33
4.3.	Safety related data	34
4.4.	Dimensions	34
4.5.	Type plate	36
5.	Installation.....	36
5.1.	Tools required.....	36
5.2.	Structure.....	37

Steripur 417

5.3.	Assembly & disassembly of the actuator.....	38
5.3.1.	Assembly.....	38
5.3.2.	Disassembly.....	39
5.4	Installation.....	40
5.4.1.	Connection butt weld ends.....	40
5.5.	Implementation.....	40
6.	Maintenance.....	41
6.1.	Wear parts.....	41
6.2.	Diaphragm change.....	42
6.2.2.	Change of two-piece diaphragms.....	43
6.3.	Cleaning.....	43
7.	Accessories.....	43
8.	Troubleshooting.....	44
9.	EC Declaration of Conformity.....	45

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Steripur 417

1. Introduction

In case you have any queries about the device, please contact our customer services stating the serial number:

SED Flow Control GmbH

Am Schafbaum 2
D-74906 Bad Rappenau
PO-Box 1306
D-74900 Bad Rappenau
Phone: +49(0)7264/921-0
Fax: +49(0)7264/921-21
E-Mail: info-sed-de@samsongroup.com
Web: sed.samsongroup.com

These operating instructions describe the technical status of the device at delivery.
It is valid for standard versions. For special versions please contact our customer services.

Any subsequent changes at the device are not considered in these operating instructions.

© 2023

The present operating instructions document is protected by copyright. Without written approval by the manufacturer it may not be translated, reproduced (mechanically or electronically) or disclosed to third parties, either in full or in parts.

The German version of these operating instructions is deciding in case of doubt, misunderstandings or translation mistakes!

Creation date: 04.02.2020
Revision: b
Printed in Germany

Steripur 417

1.1. Information about the operating instructions

Safe operation

The operating instructions contain important information for safe and correct installation of the device. Compliance with that helps preventing hazards, avoiding repair costs and downtime, and increasing reliability and operating life of the device.

Target groups

The contents of these operating instructions are intended for installation and maintenance staff.

Reading the operating instructions

All persons who work at the device must read the operating instructions in order to be familiar with correct handling and operation. The operating instructions contain important safety information! Non-compliance with such information may cause hazardous situations. The operating instructions must be read and understood.

Copyright

The present operating instructions document is protected by copyright. Without written approval by SED Flow Control GmbH it must not be reproduced, distributed, or unauthorizedly used for competitive purposes, either in full or in parts.

All rights for drawings and other documents and any power of disposal are solely with SED Flow Control GmbH; this also applies for cases of application for intellectual property rights.

Declaration of conformity

The device complies with the basic requirements of the applicable European directives. Conformity was proved.

Technical changes

We constantly strive for further development and improvement of our devices. The data contained in this edition correspond to the state of the art at the time of printing. All technical details with regard to information and figures within the operating instructions are subject to change.

Warranty

The present printed document does not contain any guarantee commitments. For that purpose, please see our general terms and conditions of delivery and payment. The prerequisite for statutory warranty is the intended use of the device in compliance with the specified operating conditions.

The warranty only applies to absence of defectiveness of the device and its components. No liability will be assumed for consequential damage occurring through failure or malfunction of the device, irrespective of their nature.

Information on the internet

Instructions and data sheets for the device are provided on the internet at:

sed.samsongroup.com

2. Safety

2.1. Safety Information

Warning!

The operating instructions contain important safety information!
Non-compliance with such information may cause hazardous situations.
The operating instructions must be read and understood.

2.2. Hazard classification

Danger!

Indicates imminent danger!
Non-compliance will result in death or severe injuries.

Warning!

Indicates a potentially hazardous situation!
Non-compliance may result in severe injuries.

Attention!

Indicates a possible hazard!
Non-compliance may result in moderate or slight injuries.

Notice!

Warning from property damage!
Non-compliance may cause damage to the device or the facility.

Indicates important additional information, tips and recommendations

Indicates important reference to information in these operating instructions and in other documentation.

Indicates an operating step that must be performed.

Steripur 417

2.3. Intended use

Diaphragm valves are made to be assembled in pipelines and are used to regulate media from high purity to the point of contaminated or even abrasive.

Use these devices only with media, the body and the sealing is consistent for.

Intended use also includes compliance with the installation and maintenance conditions prescribed by the manufacturer and the state of art.

Any other use beyond the stated purpose is regarded as not as intended. The manufacturer is not liable for any damage resulting thereof; the entire risk shall be with the user.

2.4. Misuse

- Do not supply the system with flammable or aggressive medias.
- Do not apply mechanical loads on the housing (e.g. positioning of objects on the device, using it as a step).
- Do not carry out any external modifications on the housings of devices. Do not paint any parts of the housing.
- Use the device only with components or spare parts, which are approved from SED Flow Control GmbH.
- The standard version of the Seat valve must not be used in potentially explosive areas. For this purpose please use the suitable version and use it only in approved regions!

2.5. General safety provisions

- The operating instructions by the manufacturer are binding for installation, programming, maintenance and transport.
- It is the operator's responsibility to complement the safety provisions by special instructions specific for local operating conditions, if required.
- The operating instructions and all safety-relevant rules and regulations must be kept in a safe place.
- The operating instructions and safety information must be available in full and in easily legible condition.

Accident prevention rules

In addition to the operating instructions and the binding rules and regulations for accident prevention applicable in the country of use and the place of operation, the approved technical rules for safe and professional working must be observed.

Before beginning with your work

Before beginning with your work, please obtain information about first-aid and rescue options (emergency, doctor, fire- and rescue services). Please obtain information about locations and operation of fire extinguishers and about fire alarm and firefighting options on site. Retain the unit against inadvertent operation.

During work

During operation, wear suitable work clothes.

Do not carry out activities that may impair safety.

Only operate the device in safe and functional condition.

Provisions for protection of the environment

For any work at or with the device, the provisions for environmental protection must be complied with.

Steripur 417

2.6. Residual risks

Danger!

Danger of injury through high pressure!

Pressure acting on lines and valves may cause severe injuries!

Before disconnecting any lines and valves, the pressure must be switched off and the lines must be vented.

At no time overstep the maximum operation or control pressure!

Danger through electric voltage!

Electrical voltage may cause severe injuries or death!

Before working at the facility, voltage must be switched off and secured against unauthorized activation!

Attend to valid accident prevention and safety guidelines for electrical devices!

Warning!

Protect the facility from inadvertent operation!

After shutdown the facility a controlled restart must be guaranteed!

Attention!

Danger through hot device surface while continuous operation!

Hot device surface may cause burning - or danger of fire!

Keep the device away from easy burning materials and do not touch with bare hands!

2.7. Responsibilities of the operator

Devices that are not in unobjectionable condition may cause personal injury and property damage.

It is the operator's responsibility to only operate the device in unobjectionable condition.

Hazard zones that occur between the device and any equipment on the customer's site must be secured by the operator. The operator must ensure sufficient illumination during work at the device.

2.7.1 Appointment and instruction of responsible persons

- Only employ staff who has received safety instructions.
- Responsibilities and competencies of staff with regard to installation, programming and maintenance must be clearly defined.
- Regularly monitor the staff for safe and responsible working methods and compliance with the operating instructions

2.7.2. Information obligation

It is the obligation of the operator of the device to make these operating instructions available to all persons who work with the device at any time.

Before use of the device, all persons must have read and fully understood the operating instructions.

2.8. Description of target groups

The contents of these operating instructions are intended for various target groups. The level of knowledge/information the respective target group must have is defined below.

All target groups must have read and fully understood the contents of these operating instructions.

Installation, programming and maintenance staff must

- Be 18 years of age;
- Have profound education and professional training;
- Be trained with regard to conduct in case of a fault.

3. Transport / Storage / Disposal

3.1. Delivery

3.1.1. Scope of delivery

Immediately after receipt of the delivery, please verify that the contents are not damaged and the type and scope of the delivery complies with the data stated on the delivery note.

By checking the order number, please verify that the goods received comply with the versions you ordered. In case of irregularities, please contact us immediately.

3.1.2. Notification about damage

Damage resulting from inadequate packaging or during transport must be communicated to the shipping company, the insurance, and the supplier immediately after delivery of the consignment.

3.2. Transport

Notice!

If the guidelines are not reserved, the device may be damaged!

The device must be transported in impact-resistant packaging.

During transport, the device must be protected against humidity and dirt.

The approved ambient temperature of -10 °C to $+55\text{ °C}$ must not be exceeded through the impact of heat or cold.

Steripur 417

3.3. Storage

Notice!

If the guidelines are not reserved, the device may be damaged.

Danger!

Danger of injury after reassembly!

Please check, if there are any damages and ensure, that the mounting is correct, especially loosened erection screws!

In order to keep a device that is not used over a longer period functional, a few items must be considered:

- Please store the device in its original packaging.
- The storage location must be dry and clean.
- Storage temperature must be between -10 °C and $+55\text{ °C}$
- Lock up the control air ports with protecting caps.
- Secure the device against unauthorized operation.
- Please do not store in abrasive surrounding area.

3.4. Disposal

Protect the environment!

The individual components and the packaging must be properly disposed of according to the types of materials. Please note the current rules for waste disposal!

Notice!

Please check all media contacting parts, if they are contaminated and dispose these if necessary according to the current rules of waste disposal and environmental regulations!

4. Technical Data

4.1. General technical data

Size	DN 15 – DN 50
Diaphragm Size MA	MA 25 – MA 50
Available control functions (Cf.)	Normally closed, Cf.1 & 4 Normally open, Cf.2 & 5 Double acting, Cf. 3 & 6
Orientation control air connection	Cf. 4, 5, 6, in flow direction, Cf. 1, 2, 3, 90° to flow direction
Actuator suitable for valve bodies	Two-Way bodies Welded valve configuration T-bodies Multiport bodies Tank bottom bodies
End connection valve body	Butt weld ends acc. DIN 11866 Row A (formerly DIN 11850 Reihe 1 & 2) Row B (formerly ISO 1127) Row C (formerly ASME BPE / ASTM 269) BS O.D: 4825 ; SMS 3008, JIS G 3447 Clamps, flanges and theads acc. Customer specifications or customer order
Control air connection	Threaded socket G1/4" (Standard).
Control media	Filtered, lubricated or non-lubricated air Neutral, gaseous fluids Dust- and oil content acc. DIN ISO 85731 class 4 Max. temperature 40°C
Material actuator housing	stainless steel
Material valve body	1.4435 / 316L Investment cast 1.4435 / 316L forged 1.4435 / 316L forged Fe < 0,5% (Basler Norm)
Material diaphragm	EPDM, PTFE-TM / EPDM
Installation position	Any, actuator up preferred (except tank valve)

Actuator capacity

Diaphragm Size MA	Control function 1&4:	Control function 2&5
25	0,15 NI	0,15 NI
40	0,34 NI	0,3 NI
50	0,6 NI	0,54 NI

NI: Norm litre, volume at atmospherical pressure

For Cf. 3&6 see actuator capacity of Cf.1&4 for opening and Cf.2&5 for closing.

4.2. Operating conditions

4.2.1. Operating temperatures

Warning!

Do not exceed or deceed the permitted temperatures!

The permissible temperatures are dependent on the used materials. Please pay attention to the lowest permissible temperature. For high temperatures, the maximum possible operating pressure can decrease!

Permissible media temperature for actuators

160°C dependent on application

Permissible temperature for diaphragms

Diaphragm material	Code	Media temperatur		Sterilizations temperatur
		Min.	Max.	
EPDM	18	-10 °C	90 °C	150 °C, 60 min
FPM	2	-10 °C	90 °C	Not suitable
NBR	4	-10 °C	90 °C	Not suitable
PTFE-TM / EPDM	30	-10 °C	90 °C	150 °C, 60 min
PTFE-TM / EPDM	41	-10 °C	90 °C	160 °C, 60 min, continuous operation*
PTFE-TM / EPDM	44	-10 °C	90 °C	160 °C, 60, continuous operation* min

*For continuous operation at high temperature, the valves must be regularly checked and maintained!

Permissible temperature for valve bodies

Material	Media temperatur	
	Min.	Max.
St. Steel	-10°C	160°C

4.2.2. Operating pressure

Warning!

Danger of bursting at excess pressure! Do not exceed the permitted operating pressure!

Maximal permissible pressure, unidirectional, static pressure

diaphragm	MA 25 – 50 DN 15 - 50
EPDM, FPM, NBR	10 bar
PTFE-TM / EPDM	8 bar

4.2.3. Control pressure


Warning!

Danger of bursting at excess pressure! Do not exceed the permitted operating pressure!


The control pressure has a significant impact on the lifetime of the diaphragm if you use an actuator with control function normally open or double-acting. Therefore do not use a higher control pressure as necessary.

Control function NC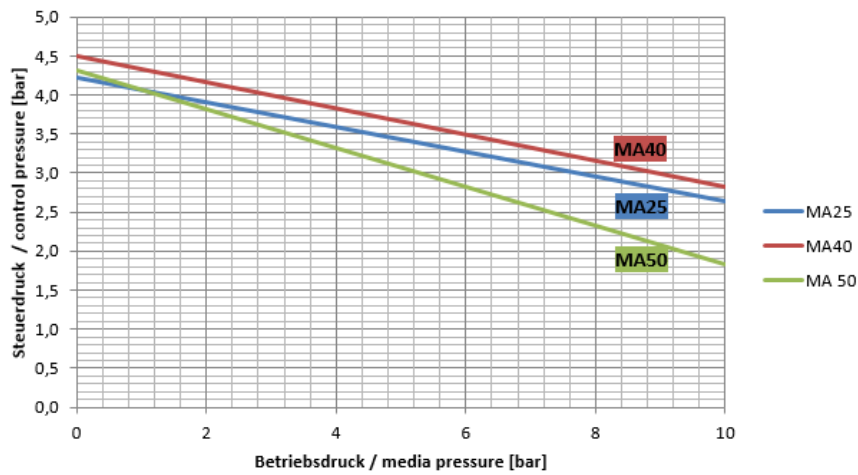

Control function DA

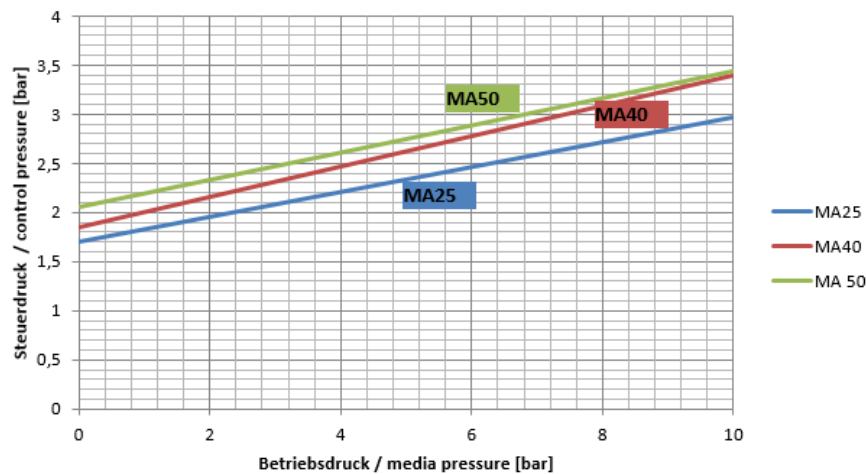Steripur 417


4.3. Safety related data

The diaphragm valve described here corresponds

- 3-A Standards Section 54-02
- TA-Luft / VDI 2440 / VDI 3479
- Pressure Equipment Directive 2014/68 / EU (declaration of conformity see attachment)

An ATEX version according to Directive 2014/34 / EU is available for use in potentially explosive areas.

4.4. Dimensions

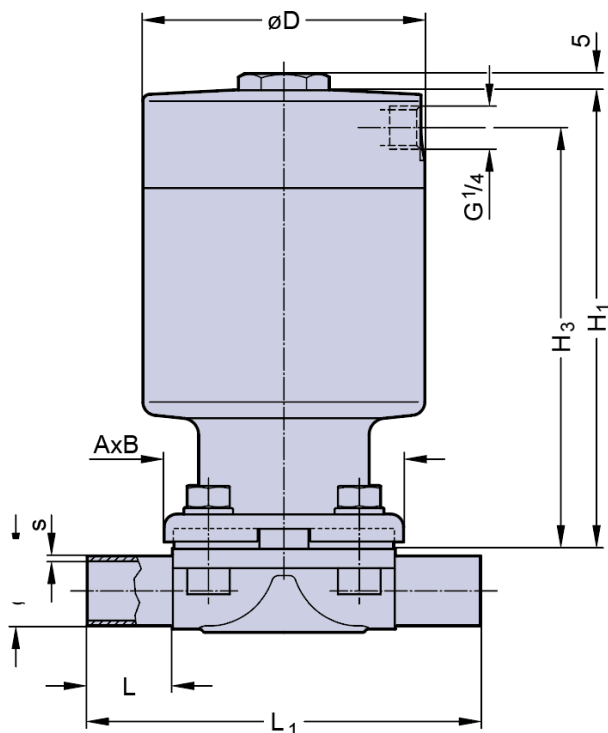DN (mm)	MA	Dimensions (mm)							Total weight ca. (kg)		Filling volume (NL)	
		L	L ₁	A x B	H ₁	H ₂	H ₃	D	Investment cast	Forged	NC	NO/DA
15-25	25	25	120	73x79	140	-	129 ^{1,2}	86	2,6	2,7	0,15	0,15
32-40	40	25	153	96x105	167	-	155 ^{1,2}	111	5,0	6,0	0,34	0,30
50	50	30	173	111x130	190	-	176 ^{1,2}	136	9,0	10,0	0,60	0,54

¹ NC, ² NO/DA

Steripur 417

Butt weld tube end standards

Butt weld tube end standard			ISO 1127 DIN 11866 Reihe B	DIN 11850		DIN Vor- zugs- reihe	ASTM 269 ASME BPE DIN 11866 Reihe C	BS O.D. 4825	SMS 3008	JIS G 3447	JIS G 3459
				Reihe 1	Reihe 2 DIN 11866 Reihe A						
Code			40	41	42	39	45 ¹	94	49	97	98 ³
DN	NPS	MA	ød x s	ød x s	ød x s	ød x s	ød x s	ød x s	ød x s	ød x s	ød x s

Valve type manually operated 205 / 206 / 290 / 297											
Valve type pneumatically operated 190 / 207 / 217											
4	-	8	-	-	-	6x1,0	-	-	-	-	-
6	-	8	-	-	8x1,0 ²	8x1,0	-	-	-	-	10,5x1,2
8	1/4	8	13,5x1,6	-	10x1,0 ²	10x1,0	6,35x0,89	-	-	-	13,8x1,65
10	3/8	8	-	12x1,0	13x1,5	-	9,53x0,89	-	-	-	-
15	1/2	8	-	-	-	-	12,7x1,65	12,7x1,2	-	-	-

Valve type manually operated 289 / 295 / 397											
Valve type pneumatically operated 188 / 195 / 317 / 392 / 394											
8	-	10	13,5x1,6	-	-	-	-	-	-	-	-
10	3/8	10	17,2x1,6	12x1,0	13x1,5	-	9,53x0,89 ³	-	-	-	17,3x1,65
15	1/2	10	21,3x1,6	18x1,0	19x1,5	18x1,5	12,7x1,65	12,7x1,2	-	-	21,7x2,1
20	3/4	10	-	-	23x1,5	22x1,5	19,05x1,65	19,05x1,2	-	-	-

Valve type manually operated 905 / 907 / 982 / 985 / 995 / 997											
Valve type pneumatically operated 385 / 395 / 402 / 407 / 417 / 495 / 592											
15	-	25	21,3x1,6	18x1,0	19x1,5	-	12,7x1,65 ³	-	-	-	21,7x2,1
20	3/4	25	26,9x1,6	22x1,0	23x1,5	-	19,05x1,65	-	-	-	27,2x2,1
25	1	25	33,7x2,0	28x1,0	29x1,5	28x1,5	25,4x1,65	-	25,0x1,2	25,4x1,2	-
32	1 1/4	40	42,4x2,0	34x1,0	35x1,5	-	31,75x1,65 ³	-	33,7x1,2	31,8x1,2 ³	-
40	1 1/2	40	48,3x2,0	40x1,0	41x1,5	-	38,1x1,65	-	38,0x1,2	38,1x1,2	-
50	2	50	60,3x2,0	52x1,0	53x1,5	-	50,8x1,65	-	51,0x1,2	50,8x1,5	-
65	2 1/2	50	-	-	-	-	63,5x1,65	-	63,5x1,6	63,5x2,0 ³	-
65	2 1/2	80	76,1x2,0	-	70x2,0	-	63,5x1,65	-	63,5x1,6	63,5x2,0 ³	-
80	3	80	88,9x2,3	-	85x2,0	-	76,2x1,65	-	76,1x1,6	76,3x2,0	-
100	4	100	114,3x2,3	-	104x2,0	-	101,6x2,11	-	101,6x2,0	101,6x2,1	-
125	5	150	139,7x2,6	-	129x2,0	-	-	-	-	-	-
150	6	150	168,3x2,6	-	154x2,0	-	152,4x2,77	-	-	-	-

Dimensions in mm; MA = diaphragm dimension / other nozzle standards upon request / preferred standards bold

¹ ASTM 269 ASME BPE tube diameter (code 45) in forged version optionally also available with tube end length according to ASME BPE (code 95); tube size 1/4" to 2 1/2" L = 1.5" (38.1 mm); tube size 3" L = 1.75" (44.45 mm); tube size 4" L = 2" (50.8 mm); tube size 6" L = 2.5" (63.5 mm).


² only valid for DIN11866

³ only available in forged design

Steripur 417

4.5. Type plate

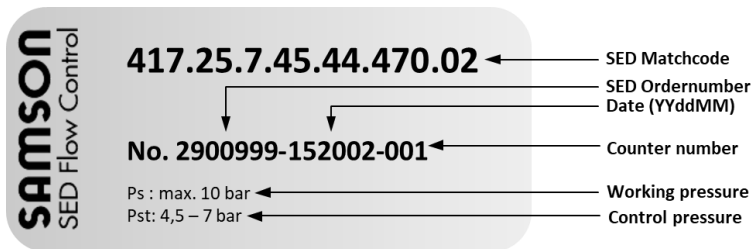The type plate is a sticker on the device and contains important information for operating. To get detailed information concerning the article codes please use the product configurator on our website, sed.samsongroup.com.

5. Installation

Danger!

Danger of injury through high pressure!

Pressure acting on lines and valves may cause severe injuries!

Before disconnecting any lines and valves, the pressure must be switched off and the lines must be vented.

Danger of injury through electric shock!

Shut off current and secure against unauthorized activation.

Warning!

Danger of injury in case of incorrect assembly!

Incorrect assembly may cause severe injuries.

Assembly and installation must only be performed by authorized expert staff.

Secure plant against unauthorized activation!

Guarantee controlled restart after shut down!

5.1. Tools required

The tools required for mounting and installation of the device are not included in the scope of delivery:


Only use functional, suitable and safe tools!

Steripur 417

5.2. Structure

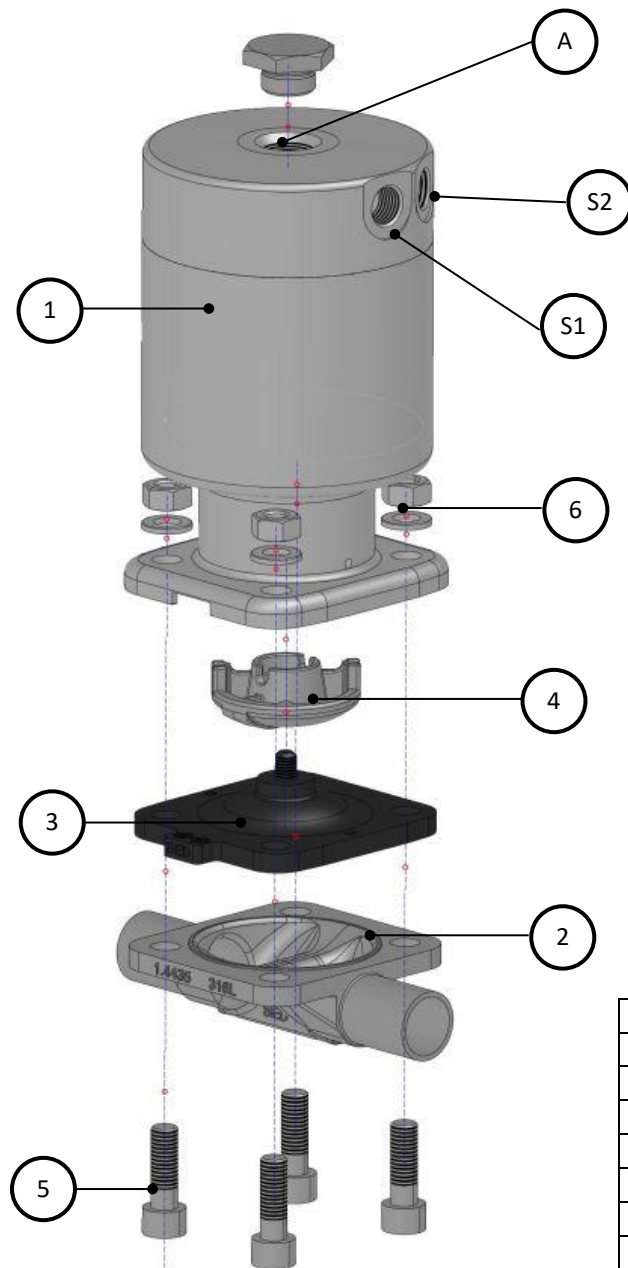1	Actuator
2	Valve body
3	Diaphragm
4	Compressor
5	Bolt
6	Washer
S1	Control air connection NC / open valve
S2	Control air connection NO / close valve
A	Assembly thread M16 for accessories

The control air connections are dependent on the control function of the actuator.

5.3. Assembly & disassembly of the actuator

Danger!

Danger of injury!

Check the device for correct assembly before implementation! Also check the tightening torque!


Only use functional, suitable and safe tools! Assembly and disassembly must only be performed by authorized expert staff.

Note!

Risk of damage !!

"In order to avoid damage inside the drive, air connection adapters to be screwed in, other screw-in adapters or pilot valves have a maximum screw-in depth of 6.5 mm. "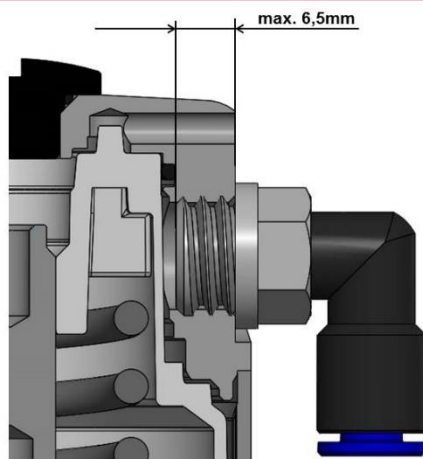

5.3.1. Assembly

For assembly of the actuator, proceed as described following:

- Open the actuator.
- Place the actuator properly on the valve body. Take care for correct orientation of the diaphragm weir and valve body weir.
- Place bolts and screw in hand-tight.
- Close pneumatic operated actuator or respectively close manual actuator by 80%.
- Fasten bolts or nuts crosswise with following tightening torque:

Steripur 417

Membrangröße <i>diaphragm size</i>	Gewindegröße <i>thread size</i>	Anzahl Schrauben <i>No. of bolts</i>	Edelstahl Ventilkörper <i>stainless steel body</i>	
			PTFE	EPDM
			Nm	Nm
MA25	M8	4	10-12	5-7
MA40	M10	4	15-17	9-11
MA50	M12	4	20-22	14-16

→ Check the valve for function and tightness
Take care for simultaneous compression of the diaphragm – about 10%, depending on diaphragm material and version. Especially PTFE-diaphragms have a setting behavior. Check the valve after the first thermal load latest and re-tighten the bolts if necessary.

5.3.2. Disassembly

Danger!

Danger of injury trough high pressure!

Pressure acting on lines and valves may cause severe injuries!

Before disconnection any lines and valves, the pressure must be switched off and the lines must be vented.

Secure plant against unauthorized activation!

Danger of injury trough media leakage (steam, hot medias, acid, lye)!

For disassembly of the actuator and diaphragm from the valve body proceed as following:

- Open the valve
- Loosen screw (5) crosswise
- Remove actuator (1) incl. diaphragm (3) from valve body (2)
- If parts are dirty, clean them. Defect parts have to be replaced by original spare parts from SED Flow Control.

For diaphragm disassembly, see chapter Diaphragm change 6.2.

5.4 Installation

Before installation, ensure that the diaphragm valve is suitable for the operating conditions. These include function, materials, operating pressure und working media in connection with operating temperature. Depending on the operating conditions and weight of the valve a suitable fixture has to be applied.

Take also care for aligning pipes.

Clean pipes before installation.

5.4.1. Connection butt weld ends

Follow the laws, directives and regulations for welding which are valid in your country.

- Disassemble the valve before welding. See chapter 5.3.2.
- Weld in valve body and cool down. If necessary, weld in in self-draining position, see chapter 5.4.1. Polish surface if needed.
- Assemble actuator and diaphragm on valve body, see chapter 5.3.1.

5.5. Implementation

Ensure correct installation before implementation. Check tightness in clamping area of actuator, diaphragm and valve body. Also check the valve for function. Control air lines have to be installed without kinks and stressless. Take care for correct connection of control air, see chapter 5.2.

6. Maintenance

Maintenance is only allowed to be carried out by trained staff. Diaphragm valves are low-maintenance parts. They have to be checked and maintained periodically. These periods are depending on working load and have to be defined according to valid rules.

6.1. Wear parts

The diaphragm is a wear part and is affected by natural wear.

The actuator is a low-maintenance part, if all regulations of this operation manual are respected.

Available spare parts are, the actuator, the media-diaphragm and the valve body.

Important Information for ordering spare parts are the article codes and order number, which you will find on the type plate of the valve.

To order spare parts, please contact us via info-sed-de@samsongroup.com or asked your local dealer.

Danger!

Danger of injury!

Bonnet is under spring pressure!

Do not open the actuator! If actuator is damaged, replace the complete actuator!

The lifetime of the diaphragm is dependent on following factors:

- Diaphragm material
- Media
- Operating temperature
- Operating pressure
- Control pressure
- Correct assembly and use

6.2. Diaphragm change

The diaphragm has to be controlled for wear after max. 100.000 (\geq MA80 50.000) switching operations. Depending on working load, the control period has to be reduced according to valid regulations and rules.

Particles in the working media can be abrasive – reduce control period accordingly.

Only use diaphragms which are suitable for your application.

Don't use higher control pressure than necessary. This helps to save the diaphragm.

Note!

Media leakage by wrong assembly!
Ensure correct assembly before startup!

Gefahr!

Danger of injury through high pressure!

Pressure acting on lines and valves may cause severe injuries!

Before disconnecting any lines and valves, the pressure must be switched off and the lines must be vented.

Protect the facility from inadvertent operation!

Danger through media leakage (steam, hot media, acid, lye)

6.2.1. Change of one-piece diaphragms

- Disassemble valve, see chapter 5.3.2.
- Bring valve in close position.
- Screw out diaphragm anticlockwise (\geq MA10).
- Insert compressor and position it correctly.

- Screw in diaphragm by hand up to the stop or respectively plug it in. Then position diaphragm by turning it back anticlockwise in a way that the wholes of the diaphragm and valve body fit together. Attention! Do not turn back more than 180°! Also take care for correct position of the diaphragm flap. If the diaphragm isn't assembled correctly, point loading can occur! This can lead to leakage and damage of the diaphragm.

- Bring valve in open position.
- Place actuator on the valve body. Take care for correct orientation of diaphragm and weir.
- Position screws and fasten hand-tight.
- Close pneumatic actuator.
- Fasten screws or nuts crosswise.

- Tightening torques see chapter 5.3.1.

- Check the valve at the pass and to the outside for tightness and function. Take care for equal compression of the diaphragm – about 10% depending on material and version. Especially PTFE diaphragms show a setting behavior. Check the valve after the first thermal load latest and re-fasten screws if necessary.

6.2.2. Change of two-piece diaphragms

- 
 - Disassemble valve, see chapter 5.3.2.
 - Bring valve in close position.
 - Screw out diaphragm anticlockwise
 - Insert compressor and position it correctly.
 - At first place the support diaphragm in correct position on the actuator and its compressor. Invert the PTFE diaphragm and screw in clockwise up to the stop. Then position diaphragm by turning it back anticlockwise in a way that the wholes of the diaphragm and valve body fit together. Attention! Do not turn back more than 180°! Also take care for correct position of the diaphragm flap.
- 
 - If the diaphragm isn't assembled correctly, point loading can occur! This can lead to leakage and damage of the diaphragm.
 - Evert PTFE diaphragm in initially position.
 - Bring valve in open position.
 - Place actuator on the valve body. Take care for correct orientation of diaphragm and weir.
 - Position screws and fasten hand-tight.
 - Close pneumatic actuator.
- 
 - Fasten screws or nuts crosswise. Tightening torques see chapter 5.3.1.
 - Check the valve at the pass and to the outside for tightness and function. Take care for equal compression of the diaphragm – about 10% depending on material and version. Especially PTFE diaphragms show a setting behavior.
- 
 - Check the valve after the first thermal load latest and re-fasten screws if necessary.

6.3. Cleaning

For external cleaning, commercially available cleaning agents can be used. Check the chemical resistance of all materials used in the valve, which could come in contact with the cleaning agent, before applying

For sterilization processes, check if the valve is suitable for the applied process.

7. Accessories

Fitting and functional accessories are available via sed.samsongroup.com or contact us.

8. Troubleshooting

Fault	Potential cause	Solution
Valve doesn't close or close fully	Control pressure too low (for NO & DA valves)	Use pressure as specified, see chapter 4.2.3.
	Media pressure too high	Do not exceed maximum control pressure, see chapter 4.2.2.
	Control media connection mixed up	Check control air connection, see chapter 5.2.
	Control media not applied	Apply control media
	Defect pilot valve	Check pilot valve for function
	Foreign matters between weir and diaphragm	Disassemble valve, remove foreign matters and check parts for damage. See Chapter 5.3.
	Defect diaphragm	Change diaphragm, see chapter 6.2.ff
	Defect actuator spring (For NC valves)	Change actuator, see chapter 5.3.ff
Valve doesn't open or open fully	Control pressure too low (for NC valves)	Use pressure as specified, see chapter 4.2.3.
	Control media connection mixed up	Check control air connection, see chapter 5.2.
	Control media not applied	Apply control media
	Defect pilot valve	Check pilot valve for function
	Wrong assembled diaphragm	Disassemble actuator and assemble diaphragm correctly. See chapter 6.2.ff
	Defect actuator spring (For NO valves)	Change actuator, see chapter 5.3.ff
Media leakage at diaphragm flange	Wrong assembled diaphragm	Disassemble actuator and assemble diaphragm correctly. See chapter 6.2.ff
	Tightening torque of screw connection actuator -diaphragm too low	Check connection of actuator and valve body, see chapter 5.3.1.
	Defect diaphragm	Change diaphragm, see chapter 6.2.ff
	Defect actuator	Change actuator, see chapter 5.3.ff
	Defect valve body	Change valve body
	Media pressure too high	Do not exceed maximum control pressure, see chapter 4.2.2.
	Operating temperature too high	Do not exceed maximum operating pressure, see chapter 4.2.1.
Media leakage at pipe connection	Defect sealing / sealing material	Change sealing / sealing material
	Improper installation	Install valve body proper
	Threaded ends loose	Re-tighten connection
Media leakage at valve body	Defect valve body	Change valve body
Control media leakage at leakage control hole	Defect stem sealing	Change actuator, see chapter 5.3.ff and check control media for pollution
Control media leakage at control air connection „open“ (for NC valves) or respectively	Defect control diaphragm or piston lip seal	Change actuator, see chapter 5.3.ff

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Bad Rappenau
HRB-Nr.: 107503
DE 411 99 172
4209/04285

Commerzbank Hellbrunn
IBAN: DE42 25 04 00 00 00 00 00 00 00 00 00
BIC: (SWIFT CODE) COBADE33HAN

SAMSON AT A GLANCE

STAFF

- Worldwide 4,500
- Europe 3,600
- Asia 600
- Americas 200
- Frankfurt am Main, Germany 1,900

MARKETS

- Chemicals and petrochemicals
- Power and energy
- District heating and cooling, building automation
- General industry
- Industrial gases
- Food and beverages
- Metallurgy and mining
- Oil and gas
- Pharmaceuticals and biotechnology
- Marine equipment
- Water and wastewater
- Pulp and paper

PRODUCTS

- Valves
- Self-operated regulators
- Actuators
- Valve accessories
- Signal converters
- Controllers and automation systems
- Sensors and thermostats
- Digital solutions

SALES SITES

- More than 50 subsidiaries in over 40 countries
- More than 200 representatives

PRODUCTION SITES

- SAMSON Germany, Frankfurt, established 1916
Total plot and production area: 150,000 m²
- SAMSON France, Lyon, established 1962
Total plot and production area: 23,400 m²
- SAMSON Turkey, Istanbul established 1984
Total plot and production area: 11,100 m²
- SAMSON USA, Baytown, TX, established 1992
Total plot and production area: 20,000 m²
- SAMSON China, Beijing, established 1998
Total plot and production area: 47,000 m²
- SAMSON India, Pune district, established 1999
Total plot and production area: 28,000 m²
- SAMSON AIR TORQUE, Bergamo, Italy
Total plot and production area: 27,000 m²
- SAMSON CERA SYSTEM, Hermsdorf, Germany
Total plot and production area: 14,700 m²
- SAMSON KTELEKTRONIK, Berlin, Germany
Total plot and production area: 1,100 m²
- SAMSON LEUSCH, Neuss, Germany
Total plot and production area: 18,400 m²
- SAMSON PFEIFFER, Kempen, Germany
Total plot and production area: 20,300 m²
- SAMSON RINGO, Zaragoza, Spain
Total plot and production area: 19,000 m²
- SAMSON SED, Bad Rappenau, Germany
Total plot and production area: 10,400 m²
- SAMSON STARLINE, Bergamo, Italy
Total plot and production area: 27,000 m²
- SAMSON VDH PRODUCTS, the Netherlands
Total plot and production area: 12,000 m²
- SAMSON VETEC, Speyer, Germany
Total plot and production area: 27,100 m²

SED Flow Control GmbH
Am Schafbaum 2 · 74906 Bad Rappenau
Telefon: +49 7264 921 0 · Telefax: +49 7264 921 21
E-Mail: info-sed-de@samsongroup.com · Internet: sed.samsongroup.com